



# MagPurix<sup>®</sup> Urine cfDNA Extraction Kit (ZP02032)

## Instructions for Use (Handbook)



Version: 1.0



48



**ZINEXTS LIFE SCIENCE CORP.**

16F., No. 93, Sec. 1, Xintai 5th Rd.,  
Xizhi Dist., New Taipei City 221416,  
Taiwan (R.O.C.)

Read and follow these Instructions for Use prior to using this product. The latest revision of this document can be found at [www.zinexts.com](http://www.zinexts.com)

# Contents

<b>Intended Use</b> .....	3
<b>Introduction</b> .....	3
<b>Kit Contents and Storage</b> .....	3
<b>Materials Required But Not Provided</b> .....	4
<b>Warnings and Precautions</b> .....	5
<b>Purification Principle</b> .....	6
<b>Before Starting</b> .....	7
<b>Preparation of sample materials</b> .....	7
<b>Procedure of MagPurix System</b> .....	8
<b>Purification Protocol - MagPurix® EVO series</b> .....	8
<b>Troubleshooting</b> .....	9
<b>Related Products</b> .....	10
<b>References</b> .....	11
<b>Symbols</b> .....	11
<b>Limited Product Warranty</b> .....	12
<b>Revision History</b> .....	12

## Intended Use

The MagPurix® Urine cfDNA Extraction Kit utilizes unique magnetic ZiBeads® technology for automated isolation and purification of circulating cell-free DNA (cfDNA) from urine samples.

The product is intended to be used by professional users, such as technicians and physicians who are trained in molecular biology techniques.

For research use only. Not for use in diagnostic procedures.

## Introduction

Product Name	MagPurix® Urine cfDNA Extraction Kit
Catalogue Number	ZP02032
Product Overview	The MagPurix® Urine cfDNA Extraction Kit is designed to extract cfDNA from Urine samples using MagPurix® EVO series automated instruments. The kit is applied with unique magnetic ZiBeads® technology, which achieves superior product quality, consistent and high product yield and reproducible results. The purified DNA is suitable for a wide range of diagnostic and research applications, including sequencing, genotyping and qPCR detection.
Applicable Instrument Model	<b>MagPurix® EVO series Instruments</b>
Display Protocol Name on The Instrument	2032 Urine cfDNA 1ml 2032 Urine cfDNA 2ml
Applicable Instrument Firmware	Check and download the latest firmware from <a href="http://www.zinexts.com">www.zinexts.com</a>
Processing Time	25-30 minutes

## Kit Contents and Storage

Shipping and Storage	The kit is shipped at room temperature. Upon receipt, store the kit at room temperature. All kit components are stable when stored properly until the expiration date shown on the kit box.	
Kit Content	The components supplied in the kit are listed below. Sufficient reagents are supplied to perform 48 purifications.	
	Contents	Amount
	<b>1</b> Reagent Cartridge	48 pcs (6x8)
	<b>2</b> Reaction Chamber	48 pcs (6x8)
	<b>3</b> Tip Holder	48 pcs (6x8)
	<b>4</b> Piercing Pin	50 pcs
	<b>5</b> Filter Tip	50 pcs
	<b>6</b> Sample Tube (7 ml)	50 pcs (25x2)

<b>7</b> Elution Tube (1.5 ml)	50 pcs
<b>S</b> Small Tip	50 pcs
Lysis Buffer 12 (500 ml)*	1 pc
Proteinase K, 10 mg/ml (10 ml)	1 pc
Zmag-S6 (1 ml) magnetic bead	2 pc
Washing Buffer 5 (100 ml)	1 pc
Barcode Sticker (EVO only)	50 pcs


Note:

\* Lysis Buffer 12 included (500 ml) is enough for sample volume up to 5 ml. Additional Lysis Buffer 12 needs to be purchased separately for sample volume >5ml.

Reagent  
Cartridge  
Contents

Each Reagent Cartridge has 10 positions with 10 sealed well.  
Positions 1-10 contain wells filled reagents for this protocol.

Reagent	Well No.
Empty	1
Empty	2
Empty	3
Empty	4
Empty	5
Empty	6
Washing Buffer 5	7
Washing Buffer 2A	8
Elution Buffer	9
Empty	10

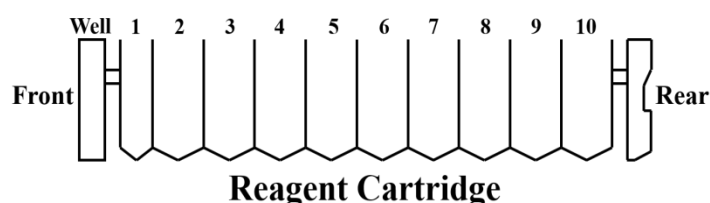

## Materials Required But Not Provided

The following general laboratory equipment and consumables are required to perform the extraction. All laboratory equipment should be installed, calibrated, operated, and maintained according to the manufacturer's recommendations. The following tables display required and special equipment along with the list of consumables.

Item
MagPurix® EVO series instrument
15 or 50 ml centrifuge tubes
Pipettes and filter tips
Phosphate-buffered saline (PBS, may be required for diluting samples)

7 ml Rack
Magnet or magnetic stand
Heat block or thermomixer with adaptor for 15/50ml centrifuge tube

## Warnings and Precautions

For *in vitro* diagnostic use only. Read all the instructions carefully before using the kit. Use of this product should be limited to trained personnel in the techniques of DNA purification. Strict compliance with the user manual is required for optimal results. Attention should be paid to expiration dates printed on the box and labels of all components. Do not use a kit after its expiration date.


When working with chemicals, always wear a suitable lab coat, disposable gloves, and protective goggles. For more information, please consult the appropriate safety data sheets (SDSs). These are available online in convenient and compact PDF format at [MSDS \(Material Safety Data Sheets\) - Downloads - www.zinexts.com](http://www.zinexts.com) where you can find, view, and print the SDSs for each kit and kit component.

Please report any serious incident occurred in relation to the device to your local representative/agent or the manufacturer, and to the competent authority of your country/state.



CAUTION: DO NOT add bleach or acidic solutions directly to the sample preparation waste.

# Purification Principle

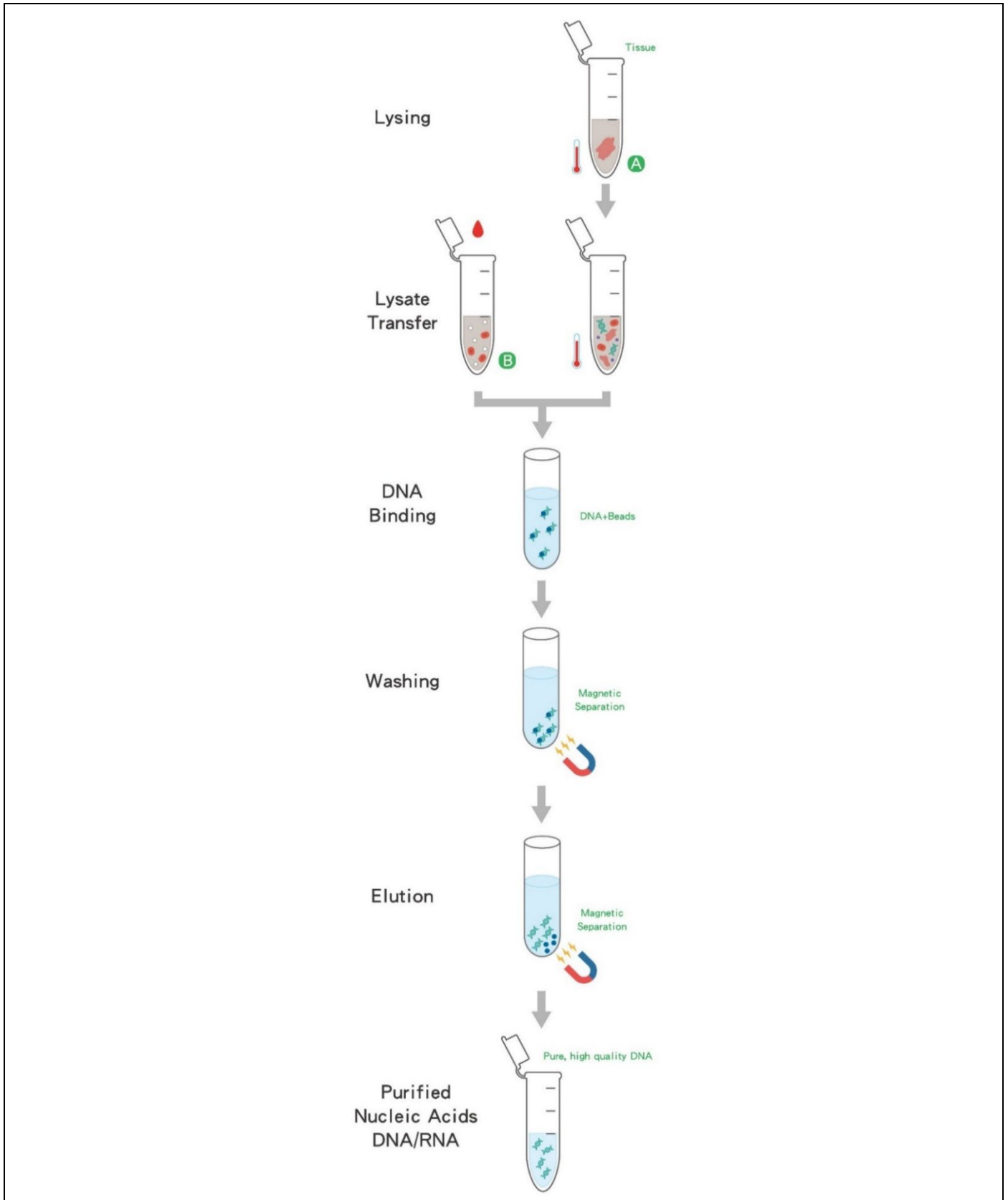

- A** Perform certain pretreatment process before extraction.
- B** Transfer sample to extraction directly.

# Before Starting

## Preparation of sample materials

The purification procedure is optimized for the use of up to 15 ml sample, such as urine, plasma, or other cell-free body fluid samples..

✂ Urine collection in the morning increases cfDNA yield as more cellular debris from the urogenital tract is shed overnight.

- a. Transfer 1-5 ml urine sample to 50 ml centrifuge tube\*
- b. Add 2 x the sample volume of Lysis Buffer 12 to sample
- c. Add 200 µl Proteinase K
- d. Add 40 µl Zmag-S6 magnetic bead
- e. Close the cap and pulse-vortex for 30 seconds.
- f. Incubate at 56°C for 30 minutes with shacking \*\*
- g. Pulse-vortex for 30 seconds
- h. Place magnet/magnetic stand on the side of the tube to hold the magnetic beads to the wall of the tube
- i. Discard supernatant while the beads held by the magnet/magnetic stand in the tube to prevent beads loss
- j. Add 1-2 ml Wash Buffer 5 to re-suspend beads
- k. Transfer mixture to 7 ml sample tube
- l. Place 7ml sample tube to 7 ml rack

### Note:

\* The Urine sample can be prepared with more than 5ml in this protocol (the maximum volume is 15ml), as long as Lysis Buffer 12 volume added in lysis step is correct (2 times the sample volume)

\*\* If the heater does not have the shacking function, perform vortex mixing every 5 minutes during incubation

The suggested starting material and elution volume range for each nucleic acid extraction.

Sample type	Starting material per sample	Elution Volume
Serum	1-5 ml (15 ml is maximum)	50-200 µl
Plasma		
Cerebrospinal Fluid (CSF)		
Urine		

# Procedure of MagPurix System

## Workflow of MagPurix operation

Place the cartridge and plastic consumables on the MagPurix instrument



Select the protocol and setup the condition



Follow onscreen message for worktable setup



Start the protocol



Collect elution product \*



UV decontamination


\* Output the bench record (option)


## Purification Protocol - MagPurix<sup>®</sup> EVO series

<b>1</b>	Turn on the Instrument	a. Turn on the power switch and wait for the screen to turn on. b. Login the instrument and enter the Home Page.
<b>2</b>	Load new Consumable(s) and Cartridge(s)	a. Open the door and remove the Sample Rack from the instrument. b. Open the Tip-Holder Lid. c. Load <b>1</b> Reagent Cartridge and all plastic disposables ( <b>2</b> Reaction Chamber, <b>3</b> Tip Holder, <b>4</b> Piercing Pins, <b>5</b> Filtered Tips and other components presented in the kit intended to use). d. Close the Tip-Holder Lid. e. Paste the barcode stickers on Elution Tubes. f. Place <b>6</b> Sample Tubes and <b>7</b> Elution Tubes into the Sample Rack.
<b>3</b>	Load the Samples	a. Transfer 7ml sample tube with mixture into the 7ml Rack. b. Put the Sample Rack back into the instrument and close the door.
<b>4</b>	Program Set up	a. Select the appropriate protocol program (for mixture volume $\leq 3$ ml select <b>2032 Urine cfDNA 1ml</b> ; for mixture volume 3-6 ml select <b>2032 Urine cfDNA 2ml</b> ) on the instrument. Press <b>NEXT</b> . b. Select the appropriate Sample Volume and Elution Volume and press <b>NEXT</b> . c. Press the number button to select the right Sample Numbers. d. Scan/Edit each primary Sample ID directly. After finished, press <b>NEXT</b> .



- e. Scan/Edit each Elution Tube ID directly. After finished, press **NEXT**.
- f. Scan Reagent Cartridge Barcode. Press **NEXT**.  
\*If the cartridge is expired, the next step cannot be performed.
- g. Follow the instructions on the screen to double-check the operating steps being completed before running the program. Press **NEXT**.

- 
- 5** Start Extraction
- a. Check “**PROGRAM CONFIRMATION**” on the screen.
  - b. Press “**START**” to start the experiment. Instrument will run the protocol program automatically until the whole process is completed.
  - c. At the end of the run (approximately 25-30 minutes), instrument alarms briefly and the screen indicates “**PROGRAM FINISH**”.
  - d. If you want to perform the same protocol, press “**RERUN**” to perform the same experiment. If you do not need to re-run the experiment, press the function button “ **HOME**” to exist the experiment mode.

- 
- 6** Collect the Elution tubes
- a. Open the instrument door.
  - b. Collect the elution tubes containing the purified nucleic acids and place on ice or in temperature lower than 8°C immediately.
  - c. The purified nucleic acids are ready for immediate use. Store the purified cfDNA at -20°C or lower temperature. However freeze-thaw cycles should be avoided, and samples should be accurately quantified once thawed prior to starting downstream analysis.
  - d. Discard the used cartridges and all plastic consumables into biohazard waste. \*Do not reuse the cartridges.
  - e. If you are not using the instrument immediately, place the Sample Rack back to the workplace, close the instrument door and press “Start” button for 2 seconds to enter sleep mode. Moreover, if the instrument will not be used in a long time, please turn off the power switch. If you are not using the instrument immediately, please put the Sample Rack back into the instrument, close the instrument door, and press the “ **POWER**” function button to enter sleep mode. If the instrument will not be used in a long time, please turn off the power switch.
- 

## Troubleshooting

\*This table is helpful for solving common problem. If you need other technical support, please contact Zinexts team (sales@zinexts.com) or your distributor.

Problem	Possible Cause	Comments and suggestions
Poor DNA quality or yield	Deterioration or contamination of reagents.	Please ensure that the kit reagents are still in the effective using period before use. Discard any kit reagent that shows discoloration or evidence of microbial contamination.
	Kit stored under non-optimal conditions	Store kit at 15-25°C at all time after arrival. If either reagent or buffer precipitate upon shipping in cold weather or during long-term storage, dissolve precipitates by gently warming and stirring the solution.

		Please do not freeze the Reagent Cartridges.
	Insufficient sample input	DNA yield depends on the sample type and the number of nucleated cells in the sample. Please proportionally adjust the total input amount of sample to increase the DNA yield.
	Too much of elution buffer was used	The elution volume can be reduced proportionally.
	The eluate of final product(s) is not enough.	Please collect issue information and provide it to your Support Representative/Technical Support as soon as possible.
Clogging issue	Too much sample material was used.	Decrease the input amount of sample material or dilute your sample.
No results in downstream analysis	No signal/The PCR was inhibited.	Using appropriate controls for analysis. Check the positive control, negative control, water (NTC) and internal control to clarify the possible causes.
Instrument malfunction/abnormal sound	Abnormal consumables: 1. Deformed Filtered Tips 2. Deformed Reaction Chamber 3. Deformed Tip Holder	Please replace the batch with normal consumables.
	Abnormal action of instrument: 1. Inaccurate correction value 2. Spare parts or components damaged	Please collect issue information (videos and pictures) and provide it to your Support Representative/Technical Support as soon as possible to calibrate or replace any other damaged or worn parts.

## Related Products










Product Name	Cat. no.
MagPurix® Blood DNA Extraction Kit 200	ZP02001
MagPurix® Blood DNA Extraction Kit 1200	ZP02002
MagPurix® Viral NA Extraction Kit	ZP02003
MagPurix® Tissue DNA Extraction Kit	ZP02004
MagPurix® Cultured Cell DNA Extraction Kit	ZP02005
MagPurix® Bacterial DNA Extraction Kit	ZP02006
MagPurix® HPV DNA Extraction Kit	ZP02007
MagPurix® TB DNA Extraction Kit	ZP02008
MagPurix® FFPE DNA Extraction Kit	ZP02009
MagPurix® Forensic DNA Extraction Kit	ZP02010
MagPurix® Viral Pathogen DNA Extraction Kit A	ZP02011
MagPurix® Viral Pathogen DNA Extraction Kit B	ZP02012
MagPurix® Viral RNA Extraction Kit	ZP02013
MagPurix® Plant DNA Extraction Kit	ZP02014
MagPurix® Total RNA Extraction Kit	ZP02015

MagPurix® Viral NA Extraction Kit LV	ZP02016
MagPurix® PLASMA CFDNA Extraction Kit	ZP02017
MagPurix® cfDNA Extraction Kit Plus	ZP02024
MagPurix® cfDNA Extraction Kit LV	ZP02025
MagPurix® Coronavirus RNA Extraction Kit	ZP02027
MagPurix® Urine cfDNA Extraction Kit	ZP02032
MagPurix® Plasma cfDNA Extraction Kit	ZP02033

## References

- Tan SC *et al.* J Biomed Biotechnol. (2009)

## Symbols

Symbol	Explanation
	For research use only
	Catalogue number
	Lot/Batch number
	Sufficient for [n] samples
	Instructions for Use
	Expiry date
	Storage temperature (15°C – 25°C)
	Manufacturer
	Caution

## Limited Product Warranty

Zinexts Life Science is committed to provide customers with high-quality products and services. Our goal is to ensure that every customer is 100% satisfied with our products and services. If you have any question or concerns, contact our Technical Support Representatives.

Zinexts Life Science guarantees the performance of all products according to the specifications stated in our product literature. The purchasers/users must determine the suitability of the product for their particular use. We reserve the right to change, alter, or modify any product to enhance its performance and design.

This warranty limits Zinexts Life Science Corporation's liability only to the cost of the product. No warranty is granted for products beyond their listed expiration date. No warranty is applicable unless all product components are stored and used in accordance with instructions.

## Revision History

Version	Date	Description
1.0	4. Nov. 2022	New edition